

EP164-EN.01

# **GDF15** [Biotinylated]: **GFRAL** Inhibitor Screening ELISA Kit

Pack Size: 96 tests

Catalog Number: EP-164

IMPORTANT: Please carefully read this manual before performing your experiment.

For Research Use Only. Not For Use In Diagnostic Or Therapeutic Procedures

FP164-FN.01

ACTO\*

**INTENDED USE** 

This kit is designed for screening of inhibitors of binding between biotinylated human GDF15 and human GFRAL.

It is intended for research use only (RUO).

PRINCIPLE OF THE ASSAY

This inhibitor screening ELISA kit is designed to facilitate the identification and characterization of new GDF15

pathway inhibitors. The assay takes advantage of our in house-developed binding of biotinylated human GDF15 and

human GFRAL in a functional ELISA assay and employs a simple colorimetric ELISA platform. Briefly, we provide

you with a human Biotinylated GDF15 protein, a human GFRAL protein, an anti-GDF15 neutralizing antibody (as

method verified Std.), and Streptavidin-HRP reagent. Your experiment will include 4 simple steps:

1) Coat the plate with human GFRAL.

2) Add your molecule of interest to the tests.

3) Add human GDF15 Biotin to bind the coated human GFRAL.

4) Add Streptavidin-HRP followed by TMB or other colorimetric HRP substrate.

Finally, the half maximal inhibitory concentration (IC50) of your compound to GDF15: GFRAL binding will be

determined by comparing OD readings among different experimental groups.

MATERIALS PROVIDED

TABLE 1. MATERIALS PROVIDED (pls modify according to COA)

2 /9



#### EP164-EN.01

Catalog	Components	Size (96 tests)	Format	Storage	
EP164-C01	High-bind Plate	1 plate	Solid	2-8	3°C
EP164-C02	Human GFRAL	10 μg	Powder	2-8°C	-70°C after
EP164-C03	Biotinylated Human GDF15	5 μg	Powder	2-8°C	reconstitution, avoid
EP164-C04	Anti-GDF15 Neutralizing Antibody	10 μg	Powder	2-8°C	freeze-thaw cycles
EP164-C05	Streptavidin-HRP	5 µg	Powder	2-8°C, avoid light	110020 11111111 0,0102
EP164-C06	Coating Buffer	12 mL	Liquid	2-8°C	
EP164-C07	20xWashing Buffer	50 mL	Liquid	2-8	3°C
EP164-C08	Blocking Buffer	50 mL	Liquid	2-8°C	
EP164-C09	Substrate Solution	12 mL	Liquid	2-8°C, avoid light	
EP164-C10	Stop Solution	7 mL	Liquid	2-8	3°C

# REAGENTS/EQUIPMENT NEEDED BUT NOT SUPPLIED

ì	Singl	e or c	iual	wavel	ength	microp	late r	eader	with	450	nm/6.	30nm	filter	,

Centrifuge;

37 °C Incubator;

Single channel or multichannel pipettes with 10 μL, 200 μL and 1000 μL precision;

 $10~\mu L$ ,  $200~\mu L$  and  $1000~\mu L$  pipette tips ;

Test Tubes;

Graduated cylinder;

Deionized or distilled water for dilution;

## **STORAGE AND VALIDITY INSTRUCTIONS**

Unopened kit should be stored at 2°C -8°C upon receiving. Find the expiration date on the outside packaging and do not use reagents past their expiration date.

The kit should be stored as TABLE 1 after the reconstitution of lyophilized materials. The shelf life is 30 days from the date of opening.

3 /9

US and Canada: Tel: +1 800-810-0816 Web: <a href="http://www.acrobiosystems.com">http://www.acrobiosystems.com</a>
Asia and Pacific: Tel: +86 400-682-2521 E-mail: <a href="mailto:order@acrobiosystems.com">order@acrobiosystems.com</a>



Note:

- a. Do not use reagents past their expiration date.
- b. Find the expiration date on the outside packaging.

## REAGENT PREPARATION

- 1. Restore all reagents and samples to room temperature (20-25°C) before use.
- 2. Reconstitute the provided lyophilized materials to stock solutions with sterile deionized water as recommended in Tab.2, Solubilize for 15 to 30 minutes at room temperature with occasional gentle mixing. Avoid vigorous shaking or vortex. The reconstituted stock solutions should be stored at -70°C. **Avoid freeze-thaw cycles**.

Note: Streptavidin-HRP stock solution should be protected from light.

TABLE 2. RECONSTITUTION METHODS FOR 96 TESTS

Catalog	Components	Amount	Stock Solution Con.	Reconstitution Buffer and Vol.
EP164-C02	Human GFRAL	10 μg	100 μg/mL	100μL, water
EP164-C03	Biotinylated Human GDF15	5 μg	100 μg/mL	50 μL, water
EP164-C04	Anti-GDF15 Neutralizing Antibody	10 μg	100 μg/mL	100 μL, water
EP164-C05	Streptavidin-HRP	5 μg	50 μg/mL	100 μL, water

# RECOMMENDED PROTOCOL

## 1. Working solution preparation

1.1 Preparation of 1×Washing Buffer:

Dilute 25 mL 20×Washing Buffer with ultrapure water/deionized water to 500 mL.

1.2 Preparation of Dilution Buffer:

10 mL Blocking Buffer (EP164-C08) add 30 mL 1×Washing Buffer.

## 2. Coating

- 1)Dilute Human GFRAL stock solution (100 μg/mL) to 0.5 μg/mL with Coating Buffer to make Human GFRAL working solution.
- 2)Add 100 μL of Human GFRAL working solution (0.5μg/mL) to each well and leave a couple of wells uncoated for No-Coating Control, seal the plate with microplate sealing film and incubate overnight (or 16 hours) at 4°C.

4 /9

US and Canada: Tel: +1 800-810-0816 Web: <a href="http://www.acrobiosystems.com">http://www.acrobiosystems.com</a>
Asia and Pacific: Tel: +86 400-682-2521 E-mail: <a href="mailto:order@acrobiosystems.com">order@acrobiosystems.com</a>

FP164-FN.01

ACTO\*

3. Washing

Remove the remaining solution by aspiration, add 300 µL of 1×Washing Buffer to each well, gently tap the plate for

1 minute, remove any remaining 1×Washing Buffer by aspirating or decanting, invert the plate and blot it against

paper towels. Repeat the washing step above for three times.

Note: For best results, the complete removal of the Human GFRAL solution is essential. The use of a manifold dispenser or an auto-

washer may be necessary.

4. Blocking

Add 300 µL Blocking Buffer to each well, seal the plate with microplate sealing film and incubate at 37 °C for 1.5

hours.

5. Washing

Repeat step 3. At the same time, you can start to prepare your samples.

6. Add Samples

1)Make serial dilution of the samples as appropriate.

2) If you intend to use the provided Anti-GDF15 Neutralizing Antibody as a reference (Std.), you may dilute the

antibody as recommended in Figure 1.

3)Add 50 µL of sample solution to each well according to our recommendation (Figure 2) or your own plate setup.

4)For No-Coating Control wells, please add 50 μL Dilution Buffer.

7.Binding

1) Dilute Biotinylated Human GDF15 stock solution (100 µg/mL) to 0.008 µg/mL with Dilution Buffer to make

Biotinylated Human GDF15 working solution.

2) For No-binding control wells, please add 50 µL Dilution Buffer.

3) For all other wells, please add 50 µL Biotinylated Human GDF15 working solution to the wells and mix the


samples by gently tapping the plate. Seal the plate with microplate sealing film and incubate at 37°C for 1 hour.

Note: The working solution should be prepared immediately before use and should not be stored.





FIG.1 PREPARATION OF 1:1 SERIAL DILUTIONS OF THE Anti-GDF15 Neutralizing Antibody

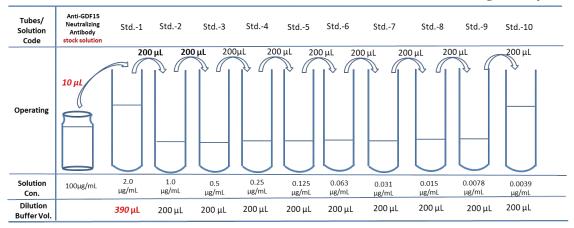

## FIG.2 PLATE LAYOUT

	1	2	3	4	5	6	7	8	9	10	11	12
Α	Std8	Std8	Std9	Std9	)				()	()	()	
В	Std7	Std7	Std10	Std10	)						(iii)	
C	Std6	Std6	Std11	Std11	)				()	()	()	
D	Std5	Std5	Std12	Std12	)	)		()	···	()	()	
E	Std4	Std4	Positive Ctrl.	Positive Ctrl.	)	)		<u></u> )	·	()	()	
F	Std3	Std3	No- binding Ctrl.	No- binding Ctrl.	X	)		····	····	()		
G	Std2	Std2	coating Ctrl.	No- coating Ctrl.	X	····)				()		
Н	Std1	Std1	coating Ctrl.	coating Ctrl.	)	)	()	()	()	()	()	

# 8. Washing

Repeat step 3.

# 9.Add Streptavidin-HRP

1)Dilute Streptavidin-HRP stock solution (50  $\mu g/mL$ ) to 0.1  $\mu g/mL$  with Dilution Buffer to make Streptavidin-HRP working solution.

6/9

**US and Canada:** Tel: +1 800-810-0816

Tel: +86 400-682-2521

Web: http://www.acrobiosystems.com

**Asia and Pacific:** 

E-mail: order@acrobiosystems.com





2)For all wells, add 100 μL Streptavidin-HRP working solution, seal the plate with microplate sealing film and incubate at 37°C for 1 hour, avoid light.

# 10.Washing

Repeat step 3.

## 11. Substrate Reaction

Add 100 µL **Substrate Solution** to each well. Seal the plate with microplate sealing film and incubate at 37°C for 20 minutes. Avoid light.

# 12.Termination

Add 50 µL Stop Solution to each well, and gently shake the plate to allow thorough mixing.

Note: the color in the wells should change from blue to yellow.

# 13.Data Recording

Read the absorbance at 450 nm using UV/Vis microplate spectrophotometer.

*Note*: Subtracting the value read at  $OD_{450nm}$  with  $OD_{630nm}$  can be used to reduce the background noise.

# **SIMPLIFIED PROTOCOL**

TABLE. 3 ASSAY PROTOCOL

Steps Code	Steps	Reagents & Instruments	Reaction Conditions	Samples	No-binding Ctrl.	No-coating Ctrl.	Positive Ctrl.
1	Working fluid preparation	N/A	N/A	N/A	N/A	N/A	N/A
2	Coating	Human GFRAL Working Solution	4°C for overnight	100 μL	100 μL	_	100 μL
3	Washing	1XWash Buffer	Wash for 3 times	300 μL	300 μL	300 μL	300 μL
4	Blocking	Blocking Buffer	37°C for 1.5 hours	300 μL	300 μL	300 μL	300 μL
5	Washing	1XWash Buffer	Wash for 3 times	300 μL	300 μL	300 μL	300 μL
	A 11 C1	Samples		50 μL	_	_	_
6	Add Samples  Dilution Buffer		_	_	50 μL	50 μL	50 μL
7	Binding	Biotinylated Human GDF15  Working Solution	Mix by gentle tapping, incubate at 37°C for 1	50 μL	_	50 μL	50 μL

7 /9

US and Canada: Tel: +1 800-810-0816 Web: <a href="http://www.acrobiosystems.com">http://www.acrobiosystems.com</a>

Asia and Pacific: Tel: +86 400-682-2521 E-mail: order@acrobiosystems.com



#### FP164-FN.01

		Dilution Buffer	hours	_	50 μL	_	_	
8	Washing	1XWash Buffer	Wash for 3 times	300 μL	300 μL	300 μL	300 μL	
9	Streptavidin-HRP	Streptavidin-HRP Working Solution	37°C for 1 hours	100 μL	100 μL	100 μL	100 μL	
10	Washing	1XWash Buffer	Wash for 3 times	300 μL	300 μL	300 μL	300 μL	
11	Substrate Reaction	Substrate Solution	37°C for 20 minutes	100 μL	100 μL	100 μL	100 μL	
12	Termination	Stop Solution	Mix by gentle tapping	50 μL	50 μL	50 μL	50 μL	
13	Data Recording	UV/Vis spectrophotometer	Measure absorbance at 450 nm, with the correction wavelength set at 630 nm					

## Note for TAB. 3:

- 1) Samples: Your samples of interest.
- 2) No-binding Ctrl.: Reaction without Biotinylated Human GDF15 added. The absorbance should be around 0.05(< 0.1) at 450 nm.
- 3) No-coating Ctrl.: Reaction without Human GFRAL coated on the wells. The absorbance should be around 0.05(< 0.1) at 450 nm.
- 4) Positive Ctrl.: Determined the max value in 450nm absorbance, when out of inhibitors.
- 5) It is recommended that all samples, controls and standards should be done in duplicates.

## **PRECAUSIONS**

- 1. This kit is for research use only and is not for use in diagnostic or therapeutic applications.
- 2. This kit should be used according to the provided instructions.
- 3. Do not mix reagents from different lots.
- 4. All reagents should be balanced to room temperature (20°C-25°C) before use.
- 5. This kit should be stored at 2°C-8°C.
- 6. Please prepare the working solution of each component according to the needs of the experiment. Except for 1x Washing Buffer, all prepared working solution is for one-time use and cannot be stored.

# **METHOD VERIFICATION**

INHIBITION OF GDF15 [Biotinylated]: GFRAL BINDING BY ANTI-GDF15 NEUTRALIZING ANTIBODY

8 /9


US and Canada: Tel: +1 800-810-0816 Web: <a href="http://www.acrobiosystems.com">http://www.acrobiosystems.com</a>

Asia and Pacific: Tel: +86 400-682-2521 E-mail: order@acrobiosystems.com



#### EP164-EN.01

Serial dilutions of Anti-GDF15 Neutralizing Antibody (Catalog # EP164-C04) (1:1 serial dilution, from 2 µg/mL to 0.000976563µg/mL) was added into GDF15 [Biotinylated]: GFRAL binding reactions. The assay was performed according to the protocol described below. Background was subtracted from data points prior to log transformation and curve fitting (QC tested).

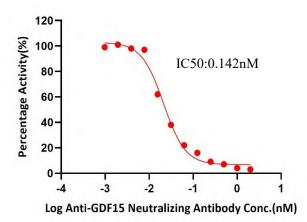

Anti-GDF15 Neutralizing Antibody conc(ug/mL)	Anti-GDF15 Neutralizing Antibody Conc(nM)	MeanAbs.(OD450)	Percentage Activity(%)
0	0.000	2.671	100%
0.001	0.007	2.656	99%
0.002	0.013	2.691	101%
0.004	0.026	2.613	98%
0.008	0.052	2.583	97%
0.016	0.104	1.658	62%
0.031	0.208	1.012	38%
0.063	0.417	0.592	22%
0.125	0.833	0.420	16%
0.250	1.667	0.249	9%
0.500	3.333	0.176	7%
1.000	6.667	0.118	4%
2.000	13.333	0.081	3%
No Coating		0.051	
No Binding		0.053	

For each experiment, a standard curve needs to be
set for each micro-plate, and the specific OD value may vary depending on different laboratories, testers, or equipments. The example data is for reference only