

Mouse Anti-RSV-F0 Antibody IgG Titer ELISA Assay Kit (Pre-Fusion glycoprotein F0)

Pack Size: 96 tests

Catalog Number: RAS-T174

IMPORTANT: Please carefully read this manual before performing your experiment.

For Research Use Only. Not For Use In Diagnostic Or Therapeutic Procedure





INTENDED USE

The kit is developed for titer measurement of Anti-RSV-F0 Antibody IgG (Pre-Fusion glycoprotein F0) in Mouse serum. It is intended for research use only (RUO).

PRINCIPLE OF THE ASSAY

Respiratory syncytial virus (RSV) is a highly contagious virus causing severe infection in infants and the elderly. Various approaches are being used to develop an effective RSV vaccine. The RSV fusion (F) subunit, particularly the cleaved trimeric pre-fusion F, is one of the most promising vaccine candidates under development. To facilitate the RSV-related research, drug trials and vaccine development, a high-throughput assay to measure IgG antibodies against the virus is in urgent need.

This assay kit is used to measure the titer of Anti-RSV-Pre-Fusion glycoprotein F0 Antibody IgG by employing an indirect ELISA. Immobilize RSV-Pre-Fusion glycoprotein F0 on the microplate. Then add the samples, incubate and wash the wells. Next add Secondary antibody HRP-Conjugated Antibody to the plate, incubate and wash the wells. Lastly load the substrate into the wells and monitor color development in proportion with the amount of antibody present. The reaction is stopped by the addition of a stop solution and the intensity of the absorbance can be measured at 450 nm and 630 nm. The OD Value reflects the amount of antibody bound.

MATERIALS PROVIDED

TABLE 1. MATERIALS PROVIDED

Catalog	Components	Size (96 tests)	Format	Storage	
				Unopened	Opened
RAS174-C01	Pre-coated RSV-Pre-Fusion glycoprotein F0 Microplate	1 plate	Solid	2-8°C	2-8°C
RAS174-C02	RSV-F0 Antibody Positive Control	100 μL	Liquid	2-8°C	2-8°C
RAS174-C03	RSV-F0 Antibody Negative Control	100 μL	Liquid	2-8°C	2-8°C
RAS174-C04	HRP-Conjugated Antibody	50 μL	Liquid	2-8°C, avoid light	2-8°C, avoid light
RAS174-C05	10×Washing Buffer	50 mL	Liquid	2-8°C	2-8°C
RAS174-C06	Dilution Buffer	50 mL	Liquid	2-8°C	2-8°C
RAS174-C07	Substrate Solution	12 mL	Liquid	2-8°C, avoid light	2-8°C, avoid light
RAS174-C08	Stop Solution	7 mL	Liquid	2-8°C	2-8°C

2/6

US and Canada:

Tel: +1 800-810-0816

Web: http://www.acrobiosystems.com

Asia and Pacific: Tel: +86 400-682-2521

E-mail: order@acrobiosystems.com





STORAGE AND VALIDITY INSTRUCTIONS

- 1. Unopened kit should be stored at 2°C-8°C upon receiving.
- 2. The opened kit should be stored per TABLE 1. The shelf life is 30 days from the date of opening.

Note: a. Do not use reagents past their expiration date.

b. Find the expiration date on the outside packaging.

MATERIALS REQUIRED BUT NOT PROVIDED

Single or dual wavelength microplate reader with 450 nm and 630 nm filter;

Centrifuge;

37 °C Incubator;

Single channel or multichannel pipettes with 10 μL, 200 μL and 1000 μL precision;

 $10 \mu L$, $200 \mu L$ and $1000 \mu L$ pipette tips;

Test Tubes:

Graduated cylinder;

Deionized or distilled water for dilution;

REAGENT PREPARATION

Bring all reagents and samples to room temperature (20°C-25°C) before use.

RECOMMENDED SAMPLE PREPARATION

1. Working fluid preparation

1.1 Preparation of 1×Washing Buffer:

Dilute 50 mL 10×Washing Buffer with ultrapure water/deionized water to 500 mL.

1.2 Preparation of Positive Control and Negative Control working fluid and pre-treatment of samples:

a. For qualitative detection of antibodies:

Dilute the samples, Positive Control and Negative Control at 1:100 with Dilution Buffer.

b. For determination of antibody titer:

It is recommended to dilute the samples, Positive Control from 1:800-1:25600 with Dilution Buffer, Negative Control from 1:100 with Dilution Buffer.

3/6

US and Canada: Tel: +1 800-810-0816 Web: http://www.acrobiosystems.com
Asia and Pacific: Tel: +86 400-682-2521 E-mail: order@acrobiosystems.com



ACTO*

2. Plate set up

Number the diluted samples corresponding to the wells of the Pre-coated RSV-Pre-Fusion glycoprotein F0 Microplate.

Each experiment requires a set of Positive Control and Negative Control working fluid.

3. Add Samples

Add 100 µL diluted samples, Positive Control and Negative Control working fluid to the corresponding wells. Add 100

μL Dilution Buffer to blank control. Seal the plate with microplate sealing film and incubate at 37°C for 1.0 h.

4. Washing

Remove the remaining solution by aspiration, add 300 µL of 1×Washing Buffer to each well, gently tap the plate for 1

min, remove any remaining 1×Washing Buffer: by aspirating or decanting, invert the plate and blot it against paper

towels. Repeat the wash step above for three times.

5. Add HRP-Conjugated Antibody

Dilute HRP-Conjugated Antibody solution at 1:1000 with Dilution Buffer to make a working solution. The prepared

working fluid should be stored away from light.

For all wells, add 100 µL HRP-Conjugated Antibody working solution. Seal the plate with microplate sealing film and

incubate at 37°C for 1.0 h, avoid light.

6. Washing

Repeat step 4.

7. Substrate Reaction

Add 100 µL Substrate Solution to each well. Seal the plate with microplate sealing film and incubate at 37°C for 20

min, avoid light.

8. Termination

Add 50 µL Stop Solution to each well, and tap the plate gently for 3 min to allow thorough mixing.

Note: the color in the wells should change from blue to yellow.

9. Data Recording

Read the absorbance at 450 nm and 630 nm using UV/Vis microplate spectrophotometer.

Note: To reduce the background noise, subtract the value read at $OD_{450 \text{ nm}}$ with the value read at $OD_{630 \text{ nm}}$.

4/6

Web: http://www.acrobiosystems.com





CUT-OFF VALUE IDENTIFICATION

Cut-off value =0.1

Normal range of Negative control (1:100): OD_{450 nm}-OD_{630 nm}< 0.1

Normal range of Positive control (1:800): OD_{450 nm}-OD_{630 nm}> 1.0

Note: The cut-off value can be determined by the end user.

INTERPRETION OF RESULTS

a. For qualitative detection of antibodies:

Positive reading: $OD_{450 \text{ nm}}$ - $OD_{630 \text{ nm}}$ of sample \geq Cut-off value means Anti-RSV-F0 Antibody IgG (Pre-Fusion glycoprotein F0) are detected.

Negative reading: $OD_{450 \text{ nm}}$ - $OD_{630 \text{ nm}}$ of sample < Cut-off value means Anti-RSV-F0 Antibody IgG (Pre-Fusion glycoprotein F0) are not detected.

b. For determination of antibody titer:

Determination of antibody titer: the positive sample was diluted with a gradient, and the antibody titer of the sample corresponds to the highest dilution factor that still yields a positive reading.

LIMITATIONS OF THE PROCEDURE

The kit cannot be used for quantitative detection.

PRECAUTIONS

- 1. This kit is for research use only and is not for use in diagnostic or therapeutic applications.
- 2. This kit should be used according to the provided instructions.
- 3. Do not mix reagents from different lots.
- 4. Bring all reagents and samples to room temperature (20°C-25°C) before use. If crystals have formed in the buffer solution, incubate until the crystals have completely dissolved. Before use, bring the solution back to room temperature.
- 5. This kit should be stored at 2°C -8°C.
- 6. Please prepare the working solution of each component according to the needs of the experiment, all prepared working solution is for one-time use and cannot be stored.

5/6

 US and Canada:
 Tel: +1 800-810-0816
 Web: http://www.acrobiosystems.com

 Asia and Pacific:
 Tel: +86 400-682-2521
 E-mail: order@acrobiosystems.com



TYPICAL DATA

Note: The Typical data is for reference only.

a. For qualitative detection of antibodies:

Value Result in units	Result	Test Result Interpretation
OD _{450 nm} - OD _{630 nm} =0.054	Negative	Anti-RSV-F0 Antibody IgG (Pre-Fusion glycoprotein F0) are not detected
OD _{450 nm} - OD _{630 nm} =0.615	Positive	Anti-RSV-F0 Antibody IgG (Pre-Fusion glycoprotein F0) are detected

b. For determination of antibody titer:

Note: Quality control data between different plates should not be mixed, and negative and positive controls should be set for each test.

Ratio of Dilution	OD _{450 nm} - OD _{630 nm} (<u>Samples</u>)	Result		
100	3.138			
200	2.977			
400	2.528			
800	1.620			
1600	0.903			
3200	0.518	The titer level of antibody is 12800		
6400	0.288			
12800	0.143			
25600	0.091			
Blank	0.016			

Web: http://www.acrobiosystems.com